

## **Stripping of Western Blot**

1. Warm up the stripping buffer to 40 °C.
2. Put the blot in stripping buffer and keep at 40 °C for 20 min.
3. Wash the blot 4 x in PBS, 20 minutes/each time.
4. Wash the blot 2 x with milk buffer, 20 minutes/each time.
5. Add 1<sup>st</sup> Ab to the blot like the regular blotting procedure.

Stripping buffer = 100ml

Tris- HCl	0.93g
SDS	2g
2-ME	200ml
pH	6.7
dH <sub>2</sub> O to	100ml